

# Freezers

Upright Laboratory & Plasma



i.Series®  
Horizon Series™

# Laboratory & Plasma Freezers

Helmer has provided high-quality laboratory equipment and refrigerated products for over 35 years. Our experience and attention to detail allow us to provide a product that offers superior uniformity, outstanding reliability, and exceptional benefits that bring value every day. In addition, our commitment to service and support ensures satisfaction throughout the life of your product. Whether storing blood products, reagents, vaccines, or medication, a Helmer product offers TrueBlue™ protection.

## Helmer Freezers Deliver

### High-Quality Craftsmanship

- Superior chamber temperature uniformity and quick recovery with heavy-duty, air-cooled refrigeration system
- Operating temperature range from -15°C to -30°C
- Bacteria-resistant powder coated interior and exterior for added security
- Sure-Seal self-closing door with magnetic closure eliminates broken door latches
- Program up to 4 daily defrost events to run during low use periods for greater efficiency
- Hospital-grade plug enhances electrical safety/serviceability

### Secure Monitoring for a Safe Storage Environment

- i.C<sup>3</sup>® with 7" (177mm) full-color touchscreen (i.Series®)
- Microprocessor temperature controller with built-in alarm/monitor (Horizon Series™)
- Key lock
- Password protected settings (i.Series)
- Chart recorder (Plasma)
- Meets AABB, ARC, FDA, EU Blood Directive 2004/33/EC, and Part 4 of the Blood Safety and Quality Regulation 2005 standards for blood plasma storage

### Designed for Convenience and Performance

- Single door models from 20 cf (572 L) to 25 cf (714 L)
- Position flush with a countertop, cabinet, or other equipment with no additional clearance required
- Recessed floor and liquid-tight drawers contain spills for easy clean-up and maintain the integrity of other products stored in the freezer
- Rechargeable battery backup (i.C<sup>3</sup>)




## i.Series® and Horizon Series™

Helmer has two distinct -30°C freezer lines. The i.Series offers many unique features not found on other freezer models. The i.C<sup>3</sup>® User Interface anchors this top-of-the-line series. It offers constant temperature monitoring and control, multiple information logs, event acknowledgement and downloading via a 7" (177mm) full-color touchscreen. (See page 5 for more information.)

The Horizon Series offers a temperature controller with a built-in alarm. It features an LED digital temperature display, programmable high and low temperature alarms, manual alarm checks, and a programmable operating range.

Both the i.Series and Horizon Series offer high quality construction and attention to detail throughout.

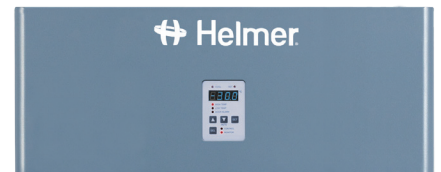

## Laboratory & Plasma Models

- i.Series and Horizon Series lines
- Capacity from 20 cf to 25 cf (572 L to 714 L)

### Upright Models

- Single door with 8 drawers or 4 shelves
- Plasma box capacity from 360 to 464 boxes
- Door-mounted, eye-level controls (i.Series)



# Upright Freezers

## Temperature Control and Monitoring

### i.Series®


- i.C<sup>3</sup>®, with 7" (177mm) full-color touchscreen monitor (described on page 5)
- Two stainless steel monitoring probes in product simulation bottles
- Temperature display in °C or °F

### Horizon Series™

- Digital microprocessor temperature controller with a built-in alarm/monitor factory set to -30°C
- Single, stainless steel monitoring probe in product simulation bottle
- Digital offset adjustment for calibration

### i.Series® and Horizon Series™

- Defrost can be programmed to actual time desired, allowing cycles to run automatically during low use periods
- Chart recorder with 4" (102mm), 7-day circular charts (Plasma)
  - Inkless, pressure-sensitive chart paper with a range of -50°C to 0°C
  - Battery backup
  - Touch button controls for chart changes and calibration

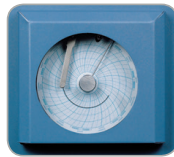

## Storage

### Laboratory Models

- Adjustable, epoxy coated wire shelves
- Stainless steel standards guard against rust and corrosion
- Side and rear guards prevent items from falling off the shelf

### Plasma Models

- 8 solid stainless steel pull out drawers with Cold-Shield™ System
- Durable, liquid-tight construction for containment of spills
- Easy-Glide stainless steel drawer slides and standards for smooth operation at low temperatures





## Refrigeration System

- Heavy-duty, air-cooled, forced-air refrigeration system with non-CFC R-404A refrigerant provides high efficiency and is environmentally friendly
- Auto condensate evaporation.
- Evaporator fan(s) shut off during door openings to maintain stable temperatures
- Programmable electric defrost

## Cabinet Construction

- Bacteria-resistant powder coated interior and exterior
- Cold-Shield™ System: Air shield with hinged door and closely spaced drawers (Plasma)
- Superior engineered frame for maximum rigidity
- Heated door frame
- Innovative self-closing Sure-Seal door system with magnetic latch
- Key lock
- Minimum of 2" (51mm) non-CFC foamed urethane insulation
- Swivel locking casters
- Access port located in top of cabinet for external monitoring probe(s)
- Rechargeable battery is environmentally friendly and long-lasting (i.C<sup>3</sup>)
- Recessed floor for easy clean-up
- Hospital-grade plug and cord

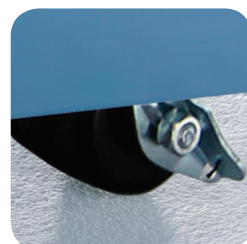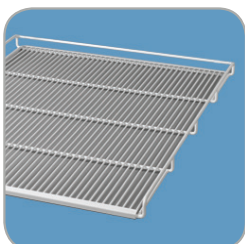



**i.C<sup>3</sup>**  
**Monitor · Acknowledge · Download**  
*Exclusively on i.Series® Models*

i.Series refrigerators are monitored with Helmer's exclusive i.C<sup>3</sup> User Interface. The i.C<sup>3</sup> provides constant temperature monitoring and control, and multiple information logs with historical data that can be downloaded and saved.


i.Act™ event acknowledgement allows users to immediately record corrective actions on-screen at the time of the event. Once downloaded, this information will be available to maintain quality control data for auditing purposes.

This icon-driven monitoring system features a door mounted 7" (177mm) full-color touchscreen ergonomically placed at eye-level on upright models and angled on undercounter models. The i.C<sup>3</sup> includes sound and visual indicators, and touch navigation. An interactive temperature graph, home screen customization, password protection, and modular design for future options are also included.


The i.D™ Integrated Electronic Access Control factory-installed option offers secure access to the unit via PIN entry on the i.C<sup>3</sup> for up to 100 users. For more information see page 8.

**i.Act™**  
**Event Acknowledgement**

Enter Event Cause, Action Taken, and Signature directly on-screen and download for audits

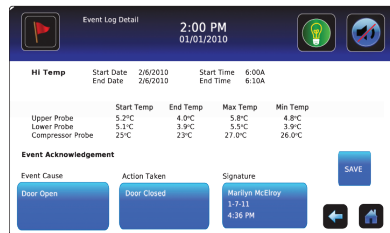

Home

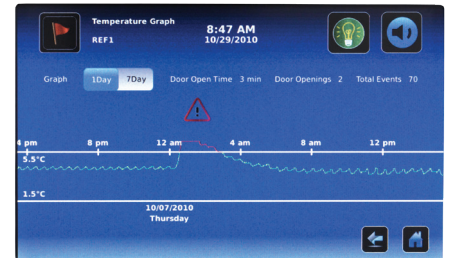



Temperature Graph



Download

**i.Series® / Horizon Series™ Monitoring**

<p><b>i.Series®</b></p> 	<p><b>Horizon Series™</b></p> 
<p><b>Exclusive i.C<sup>3</sup> User Interface</b></p> <ul style="list-style-type: none"> <li>· 7" (177mm) full-color, door-mounted touchscreen</li> <li>· Automatic high/low alarm testing (Peltier-based)</li> <li>· USB ports (2) for download/upload/upgrade</li> <li>· Optional RS232 data port</li> <li>· Dry contact alarm connection</li> <li>· Password protected settings</li> <li>· Dual alarm probes</li> <li>· Digital display of upper and lower chamber temperatures in °C and °F</li> <li>· Event log displays date, time, and temperature data</li> <li>· Compressor log displays compressor temperature performance</li> <li>· Defrost log</li> <li>· Defrost icon (snowflake) alerts user that defrost cycle is in progress</li> <li>· Interactive temperature graph</li> <li>· Min/max available through download</li> <li>· Digital calibration</li> <li>· Rechargeable battery backup</li> <li>· Date and time display</li> <li>· Alarm mute with adjustable ring-back interval</li> <li>· Adjustable alarm volume</li> <li>· Four alarm tones</li> <li>· High and low temperature alarms</li> <li>· Door ajar alarm</li> <li>· Power failure alarm</li> <li>· Condenser temperature alarm</li> <li>· RJ45 hub<sup>2</sup></li> </ul>	<p><b>Microprocessor-based controller with alarm/monitor</b></p> <ul style="list-style-type: none"> <li>· LED digital display, top bezel mounted</li> <li>· Manual alarm test</li> <li>· Dry contact alarm connection</li> <li>· Single alarm probe</li> <li>· LED display of chamber temperature in °C and °F<sup>1</sup></li> <li>· Min/max display and reset<sup>1</sup></li> <li>· Digital calibration</li> <li>· Battery backup</li> <li>· Alarm mute with 5-minute ring-back interval</li> <li>· Alarm silence with key switch</li> <li>· High and low temperature alarms</li> <li>· Door ajar alarm</li> <li>· Power failure alarm</li> </ul>

<sup>1</sup>Coming soon <sup>2</sup>For special applications

## Specifications

Model/Series	Operating Range	Volume (cu ft/L) plasma boxes	Interior Dimensions (w x h x d) (in/mm)	Exterior Dimensions* (w x h x d) (in/mm)	Shelves	Drawers	Compressor	Electrical	Net Weight
<b>Single Door</b>									
<b>iLF120</b> i.Series	-15° to -30°C	20.2 / 572	24.75 x 58.25 x 24.25 629 x 1480 x 616	29.5 x 80 x 29.5 750 x 2032 x 750	4	0	1.25 HP Air-Cooled R404A Non-CFC	208 / 230V 60Hz 4.3 FLA 230V 50Hz 3.8 FLA	443 lb / 201 kg
<b>HLF120</b> Horizon Series	-15° to -30°C	20.2 / 572	24.75 x 58.25 x 24.25 629 x 1480 x 616	29.5 x 80 x 29.5 750 x 2032 x 750	4	0	1.25 HP Air-Cooled R404A Non-CFC	208 / 230V 60Hz 4.3 FLA 230V 50Hz 3.8 FLA	440 lb / 200 kg
<b>iLF125</b> i.Series	-15° to -30°C	25.2 / 714	24.75 x 58.25 x 30.25 629 x 1480 x 769	29.5 x 80 x 35.5 750 x 2032 x 902	4	0	1.25 HP Air-Cooled R404A Non-CFC	208 / 230V 60Hz 4.3 FLA 230V 50Hz 3.8 FLA	481 lb / 219 kg
<b>HLF125</b> Horizon Series	-15° to -30°C	25.2 / 714	24.75 x 58.25 x 30.25 629 x 1480 x 769	29.5 x 80 x 35.5 750 x 2032 x 902	4	0	1.25 HP Air-Cooled R404A Non-CFC	208 / 230V 60Hz 4.3 FLA 230V 50Hz 3.8 FLA	478 lb / 217 kg
<b>iPF120</b> i.Series	-15° to -30°C	20.2 / 572 360 boxes	24.75 x 58.25 x 24.25 629 x 1480 x 616	29.5 x 80 x 29.5 750 x 2032 x 750	0	8	1.25 HP Air-Cooled R404A Non-CFC	208 / 230V 60Hz 4.3 FLA 230V 50Hz 3.8 FLA	505 lb / 230 kg
<b>HPF120</b> Horizon Series	-15° to -30°C	20.2 / 572 360 boxes	24.75 x 58.25 x 24.25 629 x 1480 x 616	29.5 x 80 x 29.5 750 x 2032 x 750	0	8	1.25 HP Air-Cooled R404A Non-CFC	208 / 230V 60Hz 4.3 FLA 230V 50Hz 3.8 FLA	502 lb / 228 kg
<b>iPF125</b> i.Series	-15° to -30°C	25.2 / 714 464 boxes	24.75 x 58.25 x 30.25 629 x 1480 x 769	29.5 x 80 x 35.5 750 x 2032 x 902	0	8	1.25 HP Air-Cooled R404A Non-CFC	208 / 230V 60Hz 4.3 FLA 230V 50Hz 3.8 FLA	557 lb / 253 kg
<b>HPF125</b> Horizon Series	-15° to -30°C	25.2 / 714 464 boxes	24.75 x 58.25 x 30.25 629 x 1480 x 769	29.5 x 80 x 35.5 750 x 2032 x 902	0	8	1.25 HP Air-Cooled R404A Non-CFC	208 / 230V 60Hz 4.3 FLA 230V 50Hz 3.8 FLA	554 lb / 252 kg

115V 60Hz NEMA 5-15 Plug  
208 / 230V 60Hz NEMA 6-15 Plug  
230V 50Hz CEE 7/7 Plug



External transformers are not utilized.

Meets IEC 61010 Standard. Meets AABB, ARC and FDA standards for blood plasma storage.

Specifications subject to change. Certified to applicable UL and CSA standards by a NRTL.

Technical Data Sheets for all Laboratory and Plasma Freezers are available at [www.helmerinc.com](http://www.helmerinc.com)

Published interior volumes are nominal; usable cu ft (liters) capacity may be offset by protrusions into cabinet.

\* Does not include door handle, electrical panel, and evaporation tray.

## Warranty

### i.Series® / Rel.i™

5 years compressor / 2 years parts / 1 year labor at Helmer

### Horizon Series™

3 years compressor / 2 years parts / 1 year labor at Helmer


Warranty applicable in US and Canada

Contact your local distributor for International warranties

## Custom Storage Options

### Roll Out Baskets

Epoxy coated wire baskets are 4" (102mm) high and adjustable in 1" (26mm) increments. Maximum load 100 lb (46 kg) per basket.

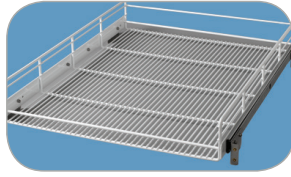


### Solid Stainless Steel Drawer with Hardware

Pull out drawers are 4" (102mm) high and adjustable in 1" (26mm) increments. Maximum load 100 lb (46 kg) per drawer.



### Full-Size Shelf with Hardware


Epoxy coated wire shelves are adjustable in 1" (26mm) increments. Maximum load 100 lb (46 kg) per shelf.

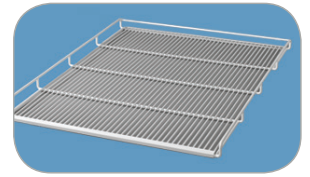

## Cabinet Options

### Temperature Chart Recorder

4" (102mm), 7-day inkless chart recorder. Includes battery backup and a supply of 52 charts. Factory installed

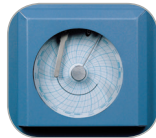

### Chart Recorder Deduction

Plasma Freezers may be ordered without chart recorder. Specify when ordering.

### Left-Hinged Single Door

Factory installed

### Stainless Steel Interior

Specify when ordering

## Accessories

### Leveling Feet

Field-installable kit. Can be installed without tipping unit.

### Storage Trays

Simplify and organize your storage needs. Each tray comes with dividers and rails so that individual storage areas can be customized. Great for storage of medication and vaccines.

### Chart Recorder Paper

4" (102mm) diameter, -50°C to 0°C, package of 52

### i.C<sup>3</sup>® Screen Protectors

Package of 3 clear screen protectors for use on i.C<sup>3</sup> User Interface

### USB Flashdrive

Used to download and upload information from/to the i.C<sup>3</sup>

### Remote Alarms (RA1 and RA4)

Extends existing alarms to remote locations within the facility. The RA1 offers monitoring of one device or zone and the RA4 provides monitoring of up to four devices or zones. Operational up to 3000 ft / 915m away (wiring not included).

### Floor and Wall Bracket Kit

Secures refrigerator to the building floor and/or wall. The unit has factory threaded mounting holes. Heavy-gauge, powder coated steel floor mounting brackets and cabinet anchor hardware are provided. Customer to provide floor and/or wall anchor hardware and onsite assembly.

### Temperature Validation

Package certifies the temperature performance of the freezer. This option must be ordered at the time the unit is purchased. Contact Helmer for additional details.


### Certificate of Calibration

Certifies that monitor was calibrated with NIST traceable thermometer before it left Helmer facility.

## Access Control

### i.Series®

#### i.D™ Integrated Electronic Access Control via i.C<sup>3</sup>® User Interface

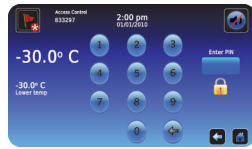

Optional Access Control offers secure access to a refrigerator or freezer. An electronic lock is integrated into

the unit and controlled by PIN entry.

- Up to 100 user ID's can be set up directly from the i.C<sup>3</sup> and programmed with an individual access PIN
- The Access Control display can be set as an alternative Home Screen.
- Key override

The Access Control log provides an audit trail of all refrigerator or freezer entries and includes the User ID, Date, Time, Duration and Method of Entry. This log can be downloaded and saved.

### Horizon Series™

#### Integrated Access Control via Keypad



The optional Horizon Access Control keypad provides fail-secure digital keypad access. A lock is integrated into the unit and unlocked by entering a user code on the integrated keypad.



- Up to 100 user codes can be entered into the keypad.
- Key override

## Also Available

### Undercounter Freezers



i.Series and Horizon Series Laboratory and Plasma Undercounter Freezers feature high-quality craftsmanship, secure monitoring, and easy access for cleaning and maintenance.

Please refer to our Undercounter Freezer brochure or [www.helmerinc.com](http://www.helmerinc.com) for more information.

