

INSTRUCTIONS MANUAL FOR:

LABS-62UCRL
FUTURA Silver Series Undercounter Refrigerator



LABRepCo, Inc
101 Witmer Road, Suite 700
Horsham, PA 19044
USA
Tel: 1-800-521-0754
Fax: 215-442-9202
www.labrepco.com
sales@labrepco.com

Table of Contents

<i>Appliance Safety</i>	3
<i>Installation Instructions</i>	4-5
<i>Appliance Use</i>	6
<i>Appliance Operation</i>	7-8
<i>Appliance Care</i>	9
<i>Temperature Selection</i>	10-11
<i>Key Lock</i>	12
<i>Alarm Instructions</i>	13-14
<i>Instructions for Built-in Applications</i>	15
<i>Instructions for Freestanding Use</i>	16
<i>Installing Panel</i>	17
<i>Before Calling for Service</i>	18
<i>Troubleshooting Guide</i>	19-20
<i>Warranty</i>	21

Appliance Safety

- Your safety and the safety of others are very important to us.
- There are many important safety messages in this manual and on your appliance. Always read and obey all safety messages.
- You can be killed or seriously injured if you do not follow these instructions.
- All safety messages will tell you what the potential hazard is, tell you how to reduce the chance of injury, and tell you what can happen if the instructions are not followed.

IMPORTANT SAFETY INSTRUCTIONS!
WARNING: To reduce the risk of fire, electric shock, or injury when using your refrigerator/freezer, follow these basic precautions:

- Plug into a grounded 3 prong outlet.
- Do not remove ground prong, nor use a 3/2 adapter
- Do not use an extension cord.
- Use care not to plug into outlets controlled by wall switches.
- Disconnect power before servicing.
- After servicing, replace all parts and panels before operating.
- Use nonflammable cleaners.
- Keep flammable materials and vapors, such as gasoline, away from appliances.
- Use two or more people to move and install appliance.
- Drawer refrigerators are provided with anti-tip brackets. Always install prior to use.

Proper Disposal of Your Old Appliance

Old appliances can represent a suffocation hazard to children

- **Remove doors from your old appliance.**
- **Leave shelves in place.**
- **If an old appliance contains refrigerants; it must be recycled by a licensed service or disposal company.**

Installation Instructions

Unpacking *Excessive Weight Hazard! * Use two or more people to move and install your appliance. Failure to do so can result in back or other injury.*

ALLOW APPLIANCE TO SIT UPRIGHT FOR 24 HOURS BEFORE TURNING ON, TO ALLOW THE COMPRESSOR LUBRICANT TO DRAIN.

1. Remove Packaging Materials
2. Remove tape and glue residue from surfaces before turning on the appliance. Rub a small amount of liquid dish soap over the adhesive with your fingers. Rinse with warm water and dry.
3. Do not use sharp instruments, rubbing alcohol, flammable fluids, or abrasive cleaners to remove tape or glue. These products can damage the surface of your appliance.
4. When Moving Your Appliance:
 - Your appliance is heavy. When moving the appliance for cleaning or service, be sure to protect the floor. Always pull the appliance straight out when moving it. Do not wiggle or “walk” the appliance when trying to move it, because floor damage can occur.
5. Clean Before Using
 - After you remove all of the packaging materials, clean the inside of your appliance before using it. See the cleaning instructions in “Appliance Care.”
6. All upright models have level legs in the front and many also have them in the rear. Level as necessary for best operation.
7. If your appliance has casters, after moving into place, lock the brakes on those casters supplied with brakes. Always unlock before moving the appliance for cleaning.

Keep flammable materials and vapors, such as gasoline, away from appliances. Failure to do so can result in explosion, or fire.

Location Requirements

- **IMPORTANT:** Do not install the appliance near an oven, radiator, or other heat source, nor in a location where the temperature will fall below 45°F (7°C). Do not place in direct sunlight.
- Check the strength of the floor before installing your appliance. It should hold the weight of a fully loaded appliance. Allow at least 1” (2.5 cm) of airspace at the top and on both sides of the appliance and 1” (2.5 cm) between the back of the appliance and the wall. If the appliance is going to be placed against a wall, if possible, leave extra space on the hinge side so the door can open wider than its width (to make it easier to remove shelves or storage bins).
- **NOTE:** This appliance may be installed side by side with another appliance. Allow for a 3” (7.5 cm) clearance at the top and sides, and a ¼” (0.64 cm) clearance between the appliances.
- The greater the airflow and ventilation you provide for your appliance—the higher its energy efficiency and performance will be.

Choosing the Right Place

- Place the appliance in a dry and regularly ventilated place. It is recommended that this appliance always be operated in an ambient temperature of at least 60°F and under 85°F. The appliance may be exposed to lower or higher temperatures but only for a short period of time and the temperature should never be lower than 45°F. Never place the appliance near heat emitting devices (e.g. cooker, radiator, water heater or similar devices) and do not expose it to direct sunlight.

Recommended Grounding Method

- A 115 Volt, 60 Hz. AC only 15- or 20- amp fused or circuit breaker, grounded electrical supply is required. It is recommended that a separate circuit serving only your appliance be provided. Use outlets that cannot be turned off by a switch or pull chain. Connect this appliance to a 3 prong power supply socket outlet (which has a ground terminal). If you only have a two prong outlet have it replaced by a qualified technician with an outlet that meets the local codes. **DO NOT USE AN EXTENSION CORD.** Required nominal voltage and frequency are indicated on the rating plate. The connection to the power supply and grounding has to be made according to current standards and regulations. The appliance resists temporary voltage tolerance from -15 to 10%.

NEVER USE INDOOR RATED APPLIANCES OUTDOORS.

Energy Saving Tips

- The more airflow you can provide around an appliance, the less energy it will use.
- If your appliance needs defrosting, keep ice build-up to less than 1/2 inch.
- Keep doors or drawers open a minimum of time, and shut securely.

Leveling Your Appliance

1. Locate the two front leveling legs taped among the packing materials. In most models these legs are already mounted into the underside of the appliance.
2. Screw these legs into the front holes on the bottom of the appliance. The back of the appliance rests on two fixed supports. (Built-in models also have rear level legs that can provide more precise leveling).
3. Adjust the front legs just enough to lift the cabinet up off of the floor.
4. Place a leveling tool on top of the appliance: first side to side, then front to back. Turn leg counterclockwise to raise a corner. Turn leg clockwise to lower a corner.
5. Repeat Step 4 until appliance is level.
6. After appliance is fully loaded, recheck with level as in Step 4.
7. If your appliance is equipped with casters, the two front casters have brakes to prevent rolling. Keep brakes locked, and unlock for moving only.

Appliance Use


- In order to ensure proper temperature, you need to permit proper airflow in the appliance. Do not block any air vents. If the vents are blocked, airflow will be prevented, and temperature and moisture problems may occur. Always position contents to allow some airflow between shelves or baskets.
- Allow the appliance to cool at least 6 to 8 hours before adding contents. Allow 24 hours for temperature to stabilize.
- NOTE: When the appliance is first started, the compressor will run constantly until the appliance is cold, It may take up to 6 or 8 hours (or longer), depending on room temperature and the number of times the appliance door is opened. After the appliance is cooled, the compressor will turn on and off as needed to keep the appliance at the proper temperature.
- The outside of your appliance may feel warm. This is normal. The appliance's design and main function is to remove heat from items and air space inside the appliance. The heat is transferred to the room air, making the outside of the appliance feel warm. *In some models the condenser that gets warm is inside the side walls, making these walls warm to the touch.*

Reversing the Door Swing (Selected models)

- If you find the direction of opening the door on your appliance inconvenient, on some models you can change it. Holes on the opposite side have already been prepared in the factory.

To Reverse Your Door

- Take off the upper hinge cover from the door and remove the screws that hold the top hinge.
- Lift the hinge straight up to free the hinge pin from the socket in the top of the door.
- Lift the door up away to free its bottom socket from the hinge pin.
- Remove the screws to the bottom hinge.
- Remove the decorative caps from opposite side and plug the holes from original side with the decorative caps.
- Move the bottom hinge to the opposite side and tighten the screws. Use any lock washers or nuts if they were installed in the original installation.
- Replace door onto the hinge pin on the bottom hinge. Place top hinge pin into door. Insert screws into the top hinge and tighten.
- If you notice the door not hanging straight—loosen the top hinges, adjust the door until it is level and tighten the screws.

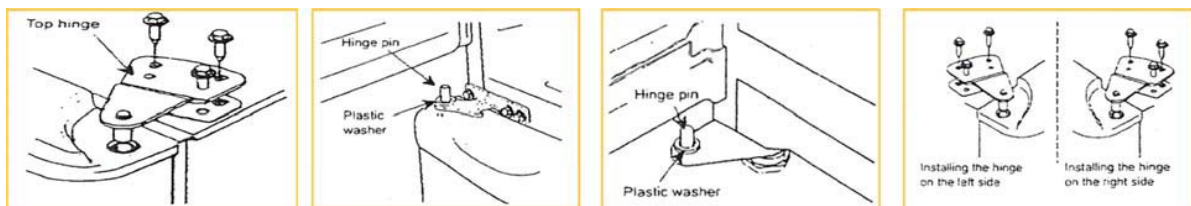

Appliance Operation

Base Grill (some models only)

To Remove and Replace the Base Grill:

1. Open the appliance door and remove the base grill by pulling down and out from the top.
 2. Replace the base grill by putting the clips in the openings in the appliance panel. Push in until the base grill snaps into place.
- An optional stainless steel grill is available. Contact your LABREPCO representative or e-mail sales@labrepc.com to find more information.

Changing the light bulb (some models only)

NOTE: Not all appliance bulbs will fit your appliance. Be sure to replace the bulb with one of the same size, shape, base size and wattage.

1. Unplug appliance or disconnect power.
2. Remove the light shield. Most models do not require tools.
3. Replace the light bulb with an appliance bulb of the same wattage.
4. Replace the light shield.
5. Plug in appliance or reconnect power.

Loss of Power

- After the power is restored, unplug the appliance for 30 minutes, to allow the refrigeration system to equalize. During this time keep the door closed whenever possible.
- **REMEMBER:** A full appliance stays cold longer than a partially filled one.

Moving Care

1. Remove all containers.
2. Unplug the appliance.
3. Clean it thoroughly. (see appliance care for cleaning instructions)
4. Tape the appliance shut and tape the electrical cord to the cabinet.

When you get to your new location, refer to the “Installation Instructions” and “Appliance Use” sections for information on installation and setting the controls.

When transporting appliances, it is recommended that you transport it upright. If it is necessary to lay it down, lay on its side—never the front or back—and be sure to check for water in the pan on top of the compressor before moving.

Defrost System, Operation and Shelf Adjustment

- These freezers are manual defrost freezers capable of operating at -25°C (-13°F). Be careful not to use sharp objects to scrape the ice. The cooling coils (part of the sealed refrigeration system) run through each fixed shelf.

- The drawers are fully removable, and you can operate the freezer without them if desired. However, the drawers contribute to energy efficiency and are best used whenever possible.
- The thermostat is located on the exterior front of the unit in the face of the tabletop. It is convenient to set it prior to installing (many customers will be satisfied with a warmer set-point, which is still colder than many household freezers).
- Some versions of this freezer have a fast-freeze button on the front of the tabletop. This function bypasses the thermostat and runs the freezer continuously (such as when a large volume of warm food is placed in the freezer). We do not recommend its use, as it wastes energy and is not necessary in most situations.
- Shelf positions for this unit are fixed in position.

Appliance Care

Cleaning Your Frost-Free Refrigerator; *Use nonflammable cleaners.*

- Frost-free means that you do not have to defrost your appliance. Frost is removed by air moving and collecting any moisture and depositing it on the cooling coil. The cooling coil periodically warms up enough to melt the frost, and the water then flows down to the drain pan (located behind the base grill) where it is evaporated.
- Complete cleaning should be done at least once a year.
 1. Unplug appliance or disconnect power.
 2. Remove all substances.
 3. Wash the inside walls with a solution of mild detergent in warm water or 2 tbs. (26 mg) baking soda to 1 qt (0.95 L) warm water.
 - Do not use abrasive or harsh cleaners such as window sprays, scouring cleansers, flammable fluids, cleaning waxes, concentrated detergents, bleaches or cleansers containing petroleum products on plastic parts, interior and door liners or gaskets. Do not use paper towels, scouring pads, or other harsh cleaning tools. These can scratch or damage materials.
 4. Rinse well and wipe dry with a clean soft cloth.
 5. Use the appropriate cleaning method for your exterior finish.
 - **Painted metal:** Wash exteriors with a clean sponge or soft cloth and a mild detergent in warm water. Do not use abrasive or harsh cleaners, or cleaners designed for stainless steel. Dry thoroughly with a soft cloth. For additional protection against damage to painted metal exteriors, apply appliance wax (or auto paste wax) with a clean, soft cloth. Do not wax plastic parts.
 6. Plug in appliance or reconnect power.
 7. Replace all containers.
 - **Stainless steel:** Clean stainless steel with special stainless steel cleaners only. Do not use bleach or any abrasive cleaners.

Cleaning Your Cycle Defrost or Manual Defrost Refrigerator

- Cycle defrost models utilize a system, where ice builds up on the rear wall and melts when the compressor is off. Manual defrost models continually build ice on inner walls.
 - **NEVER USE A METAL OBJECT TO REMOVE ICE FROM THE REAR WALL. IF NEEDED, USE A PLASTIC ICESCRAPER OR A HAIR DRYER.**
- Clean surfaces, using the same instructions as for frost-free models.
- For completely stainless steel models or models with stainless steel doors and handles; please use only cleaners especially made for stainless steel.

Temperature Selection


Your appliance may have a mechanical (dial) or digital thermostat. Instructions for both are as follows:

Mechanical Thermostat Instructions

- The refrigerator is controlled with a thermostat knob located inside the unit, towards the upper right-hand corner. Turn the knob clockwise from the **STOP (0)** position toward the higher numbers for colder temperatures.

Temperature Controls

- Higher settings on the knob (higher numbers) mean lower temperatures (colder) in all compartments of the appliance. On the coldest settings temperatures in some spots of the interior of the appliance may even drop below 32°F. Use higher thermostat settings only when required or recommended. When the ambient temperature is normal, we recommend the medium setting.
- Changes of the ambient temperature affect temperature in the interior of the appliance so choose a correct setting of the thermostat knob, depending on your ambient conditions.
- In the **STOP (0)** position the appliance does not operate (the cooling system is switched off), yet power supply is not interrupted (the light is on, when you open the refrigerator door).
- The temperature in the refrigerator and in the freezer compartment also depend on how often you open the door.
 - In all small refrigerators and freezers without interior fans, temperatures are coldest near the bottom, and occasionally near the rear. Place contents accordingly. It is best not to operate these units mostly empty. If temperatures are erratic, try putting some water bottles or freezer packs inside to create a load for the compressor.

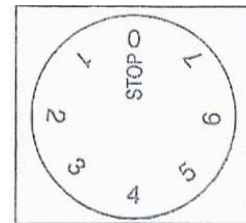

Digital Thermostat Instructions

- If you find that it is necessary to change the temperature setting for your digital thermostat controlled appliance, please refer to the following instructions. The front panel of the thermostat is illustrated here:

To see the *SETPOINT*:

- Push and immediately release the *SET* key -the display will show the *SET* point value. Wait 5 seconds and the probe value will display again.

To change the *SETPOINT*:

- Push and hold the *SET* key for more than 3 seconds to change the *set point* value; The value of the *set point* will be displayed, and the “snowflake” LED will start blinking.

To change the *SET* value:

- Push the up or down arrows to the desired setting; the new setting will be memorized after 15 seconds, provided no other keys are pushed.
- Note: The minimum and maximum range of your thermostat is preset and locked by us; so it cannot easily be accidentally moved between freezer and refrigerator temperatures. To reset this range lock, either call us or find information on our website under troubleshooting.

Key Lock (optional on most models)

Front Lock


- Your lock has a feature called automatic key eject. To lock or unlock the appliance, you must push and hold the key in the lock while turning. The lock is spring loaded, so the key cannot be left in the lock.
- If the key is lost, contact your LABREPCO representative or e-mail sales@labrepc.com to purchase a replacement. Always give the model and serial number of your appliance when requesting a replacement key.

***NEVER KEEP KEYS WITHIN THE REACH OF CHILDREN**

Alarm Instructions (alarms are optional on most models)

Setting the High/Low Alarm

1. Press the MODE button until unit is in the ALARM DISPLAY MODE indicated by the display of current temperature and HI and LO alarm settings.
2. To set the HIGH temperature alarm limit press the HI button, to advance the display in 1° increments, until the desired high temperature is displayed. Press and hold the HI button to display in 5° increments.
3. To set the LOW temperature alarm limit press the LO button, to advance the display in 1° increments, until the desired low temperature is displayed. Press and hold the LO button to rapidly advance the display in 5° increments.

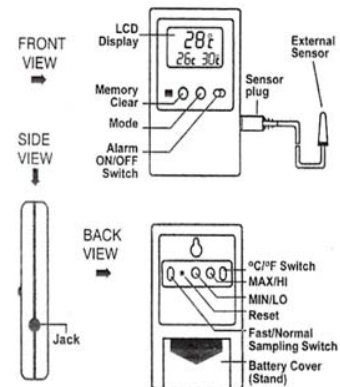

Alarm

- The unit will sound the alarm in both NORMAL DISPLAY MODE and ALARM DISPLAY MODE. Slide the ALARM ON/OFF switch to ON. When the current temperature display rises above the HI set point, or falls below the LO set point, the alarm will sound for 1 minute and issue a 3 second repeater beep every minute thereafter, for up to 12 hours. The alarm will continue to sound even if the current temperature display returns to an “in range” condition.
- To temporarily silence the alarm press either the HI or LO button on the back of the unit. The unit is still active and the alarm will sound again if the temperature goes out of range. To permanently silence the alarm, slide the ALARM ON/OFF switch to OFF position.

Operation Modes

- These alarms have two operating modes: NORMAL AND DISPLAY
 - DISPLAY MODE and ALARM DISPLAY MODE. To change from one mode to another press the MODE button once.
 - NORMAL DISPLAY MODE - is indicated by the display of current temperature and the MIN (minimum) and MAX (maximum) temperatures. These temperatures relate to external (probe) temperature if the probe is plugged in, and to ambient temperature if the sensor probe is removed.
 - ALARM DISPLAY MODE - is indicated by the display of current temperature and the HI and LO alarm settings. These temperatures relate to external (probe) temperature if the probe is plugged in, and to ambient temperature if the sensor probe is removed.

Resetting the Unit

- Use a pointed object to push the RESET button on the back of the unit. It is necessary to reset the unit whenever a change is made to °C/°F temperature, FAST/NORMAL sampling rate, or the probe sensor is plugged in or removed.

- **NOTE:** HI/LO temperature alarm settings and MIN/MAX memories will be cleared after resetting the unit.

Selecting the Sample Rate

- To select normal sampling rate (60 seconds) or fast sampling rate (10 seconds) slide the switch on back of unit to NORMAL or FAST
- **NOTE:** The unit must be reset after changing the sampling rate. Using the NORMAL sampling rate will extend battery life.

Resetting Min/Max Memories

There are two ways to reset the memories:


1. To reset both minimum and maximum memories - press MEMORY CLEAR button.
2. To reset individual minimum and maximum memories - press either the MIN or MAX button (on back of unit) once to reset the respective memory. The other memory will remain unaffected.

NOTE: You must be in Normal Display Mode to reset the memories.

Instructions for Built-In Applications


Note: The directions below are for installation under a cabinet. If you will be using this product in a freestanding application, begin with step A1 in lower half of page.

1. Remove door from chassis by unscrewing 3 screws in hinge. Keep screws, discard hinges.
2. Remove top and bottom hinge bushings.
3. Remove frame from door.
4. Attach customer supplied handle to door panel using 3 screws.
5. Screw door panel to frame using 14 screws.
6. Slide frame over refrigerator door and attach with 1 screw on each side on center of door.
7. Attach top brace to frame with 1 screw on each side. Attach bottom brace in the same manner.
8. Drill a hole through the top and bottom panel using the braces as a guide with a 1/8" drill bit. Install screws to finish attaching the top and bottom braces securely to the frame.
9. Drill a hole into the door panel from the top and bottom using the braces as a guide with 1/4" drill bit. Do not exceed 5/8" deep. This is the hole that the door will pivot on.
10. Remount the door using the supplied top and bottom hinges.

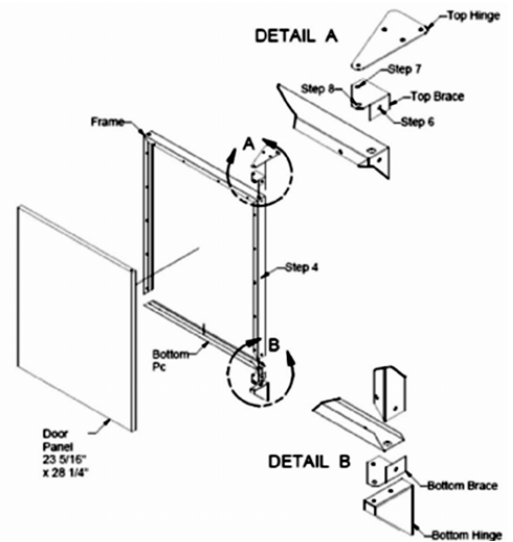

Instructions for Freestanding Use

- A1. Remove door from chassis by unscrewing 3 screws in top hinge and 2 screws in bottom hinge. Keep screws, discard hinges.
- A2. Remove top and bottom hinge bushings.
- A3. Remove frame from door.
- A4. Attach customer supplied handle to door panel using countersunk screws.
- A5. Screw door panel to frame using 14 screws.
- A6. Slide frame over refrigerator door and attach with 3 screw on each side
- A7. Install top and bottom hinge bushings.
- A8. Remount the door using the supplied top and bottom hinges.

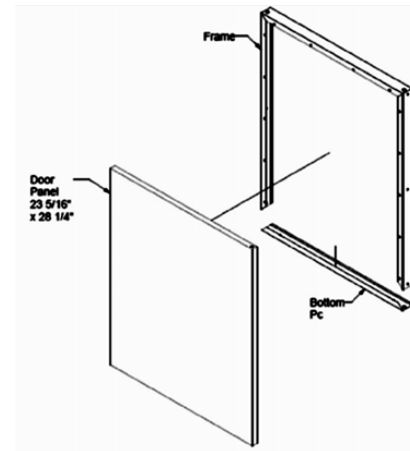

INTEGRATED FRAMES ARE OPTIONAL ON ALL MODELS.

Installing a Panel

(optional on most counter height models)

Your refrigerator or freezer is fully finished and does not need a front panel. However, if you have ordered a model with our optional frame, follow the directions below to install your panel.

- 1) Prepare your panel. Panel size is 23 1/8" W x 28 1/8" H by 1/4" thick at the edges)
- 2) Carefully lay the refrigerator (it must be empty) on its back.
- 3) Unscrew the small piece of frame from the bottom. For a thin metal panel use a backing piece of cardboard
- 4) Slide panel into frame and replace the bottom frame and reattach the screws.
- 5) Lift refrigerator upright carefully.

Before Calling for Service

- Before calling for assistance or service, please check “Troubleshooting.” It may save you the cost of a service call. If you still need help, follow the instructions below.
- When calling, please know the purchase date and the complete model and serial number of your appliance. This information will help us to better respond to your request.
 - 1) Check the plug, fuse and/or circuit breaker.
 - 2) Test the outlet with a lamp
 - 3) Make certain that the outlet used is not controlled by a wall switch. This is more common in outdoor kitchens.
 - 4) Be sure the control is not set to **0 (OFF)**.
 - 5) On outdoor units, check that the GFCI switch on the line cord is on. If necessary press the reset button.
 - 6) Read the manual, especially the items under troubleshooting. You may find the answer to your question.

If you don't find the answer to your question in this manual, contact your LABREPCO representative or e-mail sales@labrepc.com.

If You Need Replacement Parts

- If you need to order replacement parts, contact your LABREPCO representative or e-mail sales@labrepc.com.

Troubleshooting

Try the solutions suggested here first to avoid the cost of an unnecessary service call...

Your refrigerator or freezer will not operate:

- Has a household fuse blown or circuit breaker tripped, or is the outlet not receiving power or switch controlled?
 - Replace the fuse or reset the circuit breaker. Check the outlet with a small appliance you know works.
- Is the refrigerator or freezer defrosting (on some models)?
 - Recheck to see whether the appliance is operating in a few hours. It may regularly run an automatic defrost cycle.

There is water in the defrost drain pan (on some models):

- Is the appliance defrosting?
 - The water will evaporate. It is normal for water to drip into the defrost pan.
- Is it more humid than normal?
 - When it is humid, expect that the water in the defrost pan will take longer to evaporate.

The motor seems to run too much:

- Is there excessive frost or a package preventing the door from closing?
 - Defrost and clean the appliance, or move the package so the door closes properly.
- Is the room temperature hotter than normal?
 - The motor will run longer under warm conditions. At normal room temperatures, your motor will run about 40% to 80% of the time. Under warmer conditions, it will run even more.
- Has a large amount of food just been added to the appliance?
 - Adding a large amount of food warms the refrigerator or freezer. The motor normally will run longer to cool the appliance back down.
- Is the door closed completely?
 - Push the door firmly shut. If it will not shut all the way, see “The door will not close completely” later in this section.
- Are the condenser coils dirty (on some models)?
 - This prevents air transfer and makes the motor work harder. Clean the condenser coils.
- Is the door gasket sealed all the way around?
 - Contact a qualified technician.

NOTE: If the problem is not due to any of the above, remember that the compressor on your new appliance will operate longer than your old one due to its high-efficiency motor.

The interior temperature is too warm:

- Are the air vents blocked?
 - This prevents cold air movement. Move objects away from the air vents. See “Ensuring Proper Air Circulation” for air vent locations.
- Is the door opened often?
 - Be aware that the refrigerator or freezer will warm when this occurs. In order to keep the appliance cool, try to get everything you need out of it at once, keep food organized so it is easy to find, and close the door as soon as the food is removed.
- Has a large amount of food just been added?
 - Adding a large amount of food warms the refrigerator or freezer. It can take several hours to return to the normal temperature.

The door will not close completely or is difficult to open:

- Are food packages blocking the door?
 - Rearrange containers so that they fit more tightly and take up less space.
- Are the shelves and basket out of position (on some models)?
 - Put the shelves and basket back into their correct positions.

Specifications are subject to change without notice.

Warranty

Call customer service toll free in the U.S.A. @ 1-800-521-0754; Monday to Friday 8:30 am. - 5:00 p.m. (ET).

ONE-YEAR LIMITED WARRANTY

Within the 48 contiguous United States, for one year from the date of purchase, when this appliance is operated and maintained according to instructions attached to or furnished with the product, warrantor will pay for factory-specified parts and repair labor to correct defects in materials or workmanship. Service must be provided by a designated service company. Outside the 48 states, all parts are warranted for one year from manufacturing defects. Plastic parts, shelves and cabinets are warranted to be manufactured to commercially acceptable standards, and are not covered from damage during handling or breakage.

5-YEAR COMPRESSOR WARRANTY

- The compressor is covered for 5 years.
- Replacement does not include labor.

ITEMS WARRANTOR WILL NOT PAY FOR:

- Service calls to correct the installation of your appliance, to instruct you how to use your appliance, to replace or repair fuses or to correct wiring or plumbing.
- Service calls to repair or replace appliance light bulbs or broken glass shelves. Consumable parts (such as filters) are excluded from warranty coverage.
- Damage resulting from accident, alteration, misuse, abuse, fire, flood, acts of God, improper installation, installation not in accordance with electrical or plumbing codes, or use of products not approved by warrantor.
- Replacement parts or repair labor costs for units operated outside the United States.
- Repairs to parts or systems resulting from unauthorized modifications made to the appliance.
- Expenses for travel and transportation for product service in remote locations.
- The removal and reinstallation of your appliance if it is installed in an inaccessible location or is not installed in accordance with published installation instructions.

DISCLAIMER OF IMPLIED WARRANTIES; LIMITATION OF REMEDIES

CUSTOMER'S SOLE AND EXCLUSIVE REMEDY UNDER THIS LIMITED WARRANTY SHALL BE PRODUCT REPAIR AS PROVIDED HEREIN. IMPLIED WARRANTIES, INCLUDING WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE, ARE LIMITED TO ONE YEAR. WARRANTOR SHALL NOT BE LIABLE FOR INCIDENTAL OR CONSEQUENTIAL DAMAGES. SOME STATES DO NOT ALLOW THE EXCLUSION OR LIMITATION OF INCIDENTAL OR CONSEQUENTIAL DAMAGES, OR LIMITATIONS ON THE DURATION OF IMPLIED WARRANTIES OF MERCHANTABILITY OR FITNESS, SO THESE EXCLUSIONS OR LIMITATIONS MAY NOT APPLY TO YOU. THIS WARRANTY GIVES YOU SPECIFIC LEGAL RIGHTS AND YOU MAY ALSO HAVE OTHER RIGHTS, WHICH VARY FROM STATE TO STATE.

LABRepCo, Inc
101 Witmer Road, Suite 700
Horsham, PA 19044 USA
Tel: 1-800-521-0754
Fax: 215-442-9202
www.labrepco.com

LABRepCo
One Company • Many Solutions
For parts and accessory ordering,
troubleshooting and helpful hints, visit: