

STAINLESS STEEL ANGULAR RACKS WITH COLLECTION PANS FOR KEYHOLE SUSPENDED PRIMATE CAGES

Series O Racks fit Series O Cages



100123-00 Cages (shown)

104077-00 Rack (shown)

This Stainless Steel Angular Rack with Collection Pans for Keyhole Suspended Primate Cages is fabricated from die-formed angles and channels. The angle uprights' lower ends have caster couplers. The angle uprights, angle, and channel crossmembers are butted at all joints and heli-arc welded for strength. Each cage is held steadily in place with three keyhole interlocking studs.


The rack has a one-piece deflector, consisting of back and sides, that directs all wastes into the removable tapered collection pans. The pans are supported by full-length runners that are heli-arc welded to the rack's crossmembers. In addition, the pans are locked in place with a pivot type latch.

The rack has a glass beaded satin finish, removing any discoloration.

Racks are equipped with chrome plated swivel casters with 5" polyolefin wheels, two with wheel brakes. Swivel and axle bearings have seals and zerk grease fittings.

OPTIONS

Stainless Steel watering system
CPVC watering system
Stainless Steel casters
Floor Levelers

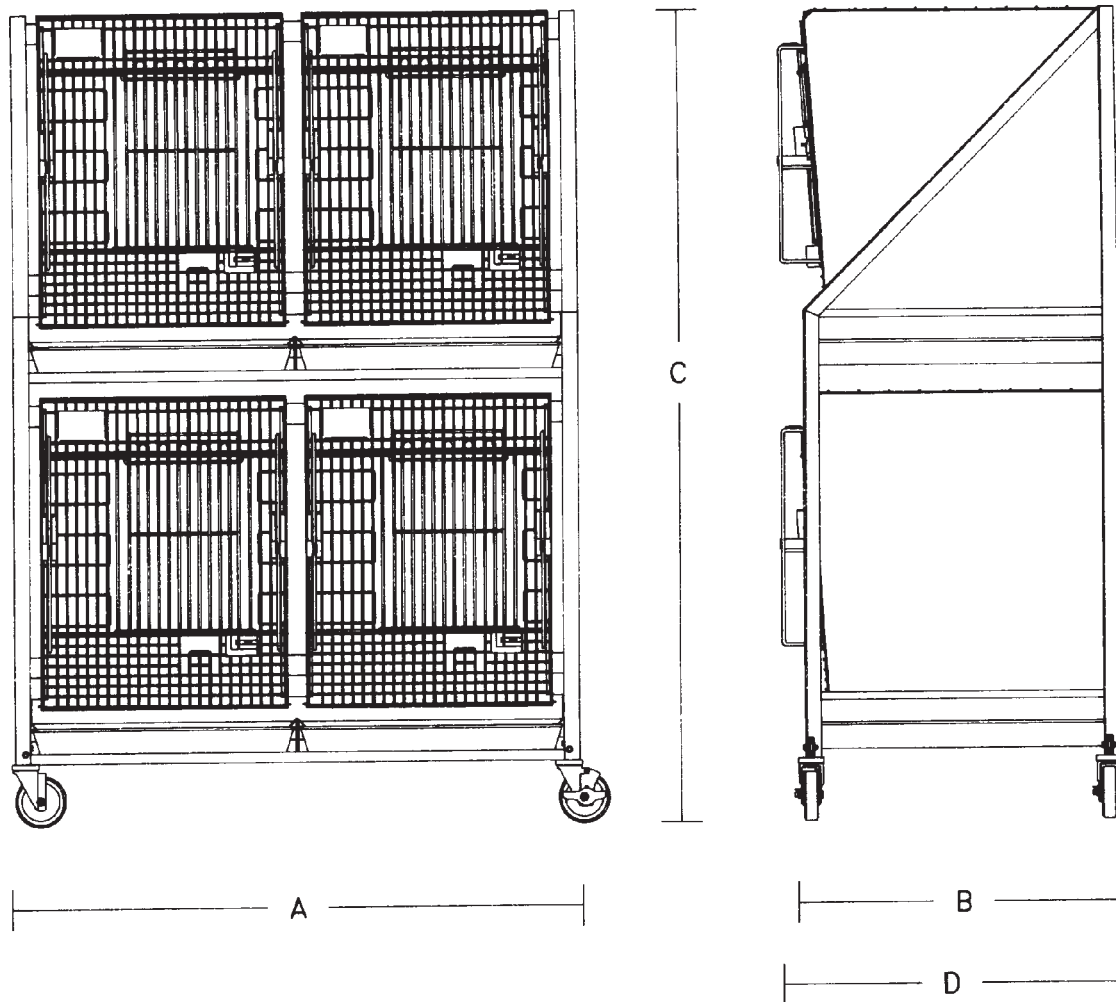

**STAINLESS STEEL ANGULAR RACKS WITH COLLECTION PANS
FOR KEYHOLE SUSPENDED PRIMATE CAGES**

PRODUCT NUMBER	CAGES	RACK CAPACITY		RACK DIMENSIONS				& 1 SLIDING DOOR (A)	& 2 SLIDING DOORS (A)
		CAGES WIDE	TIERS HIGH	WITHOUT CAGE	WITH CAGE	C	D		
For 1.6 square foot cages									
104019-00	4	2	2	36.000"	24.938"	66.500"	26.250"		
104020-00	6	3	2	52.000"	24.938"	66.500"	26.250"		
104021-00	8	4	2	68.000"	24.938"	66.500"	26.250"		
For 3 square foot cages									
104045-00	2	1	2	26.000"	32.125"	79.500"	33.438"		
104049-00	4	2	2	48.000"	32.125"	79.500"	33.438"		
104050-00	6	3	2	70.000"	32.125"	79.500"	33.438"		
For 4.3 square foot cages									
104073-00	2	1	2	30.000"	32.125"	79.500"	33.438"		
104077-00	4	2	2	56.000"	32.125"	79.500"	33.438"		
104078-00	6	3	2	82.000"	32.125"	79.500"	33.438"		
For 6 square foot cages									
104129-00	2	1	2	40.000"	32.125"	83.500"	33.438"		
104133-00	4	2	2	76.000"	32.125"	83.500"	33.438"		

ADDITIONAL EQUIPMENT

Stainless steel tapered collection pans

601673 for 1.6 sq ft cage

602995 for 3 sq ft cage

608897 for 4.3 sq ft cage

608069 for 6 sq ft cage

Racks with other cage capacities available

STAINLESS STEEL ANGULAR FLUSH RACKS WITH DOWNSPOUTS FOR KEYHOLE SUSPENDED PRIMATE CAGES

Series O Racks support Series O Cages



100133-00 Cages (shown)

104071-00 Rack (shown)

This Stainless Steel Angular Flush Rack with Downspouts for Keyhole Suspended Primate Cages is fabricated from die-formed angles and channels. The angle uprights' lower ends have caster couplers. The angle uprights, angle crossmembers, and channel crossmembers are butted at all joints and welded for strength. Each cage is held securely in place with three keyhole interlocking studs.


Under each tier of cages is a stainless steel welded trough pitched towards the downspout. The trough's ends and back are formed-up above the bottoms of the suspended cages. In addition, the trough's front edge is formed-up to create a lip.

The rack has a glass beaded satin finish, removing any discoloration.

The rack is equipped with chrome plated swivel casters with 5" polyolefin wheels, two with wheel brakes. Swivel and axle bearings have seals and zerk grease fittings

OPTIONS

Stainless Steel watering system
CPVC watering system
Stainless Steel casters
Floor levelers

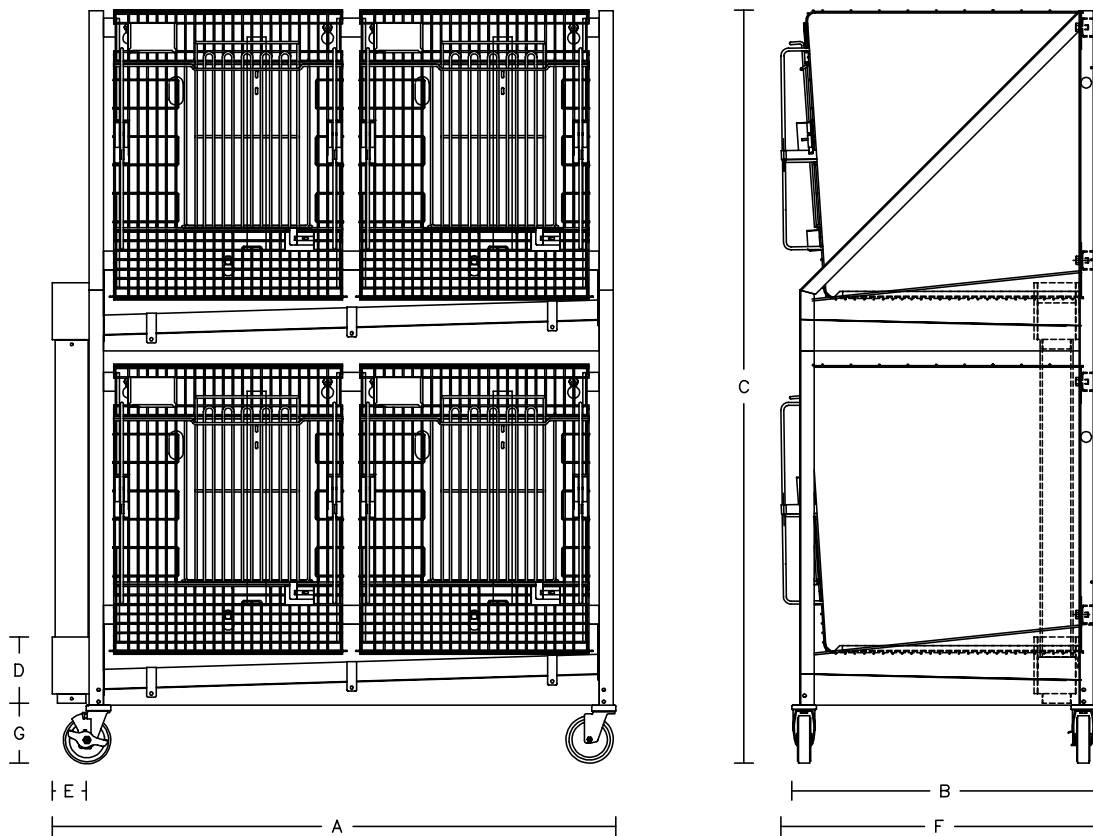


**STAINLESS STEEL ANGULAR FLUSH RACKS WITH END DOWNSPOUTS
FOR KEYHOLE SUSPENDED PRIMATE CAGES***

PRODUCT NUMBER	SINGLE WIDTH CAGES	RACK CAPACITY		RACK DIMENSIONS							
		CAGES WIDE	TIERS HIGH	A	B	WITHOUT CAGE C	D	E	WITH CAGE F	G	
For 1.6 square foot cages											
104015-00	4	2	2	40.938"	24.938"	66.718"	6.250"	5.250"	26.250"	6.375"	
104016-00	6	3	2	56.938"	24.938"	66.718"	6.250"	5.250"	26.250"	6.375"	
104017-00	8	4	2	72.938"	24.938"	66.718"	6.250"	5.250"	26.250"	6.375"	
For 3 square foot cages											
104038-00	2	1	2	30.938"	32.125"	79.500"	6.250"	5.250"	33.438"	6.375"	
104042-00	4	2	2	52.938"	32.125"	79.500"	6.250"	5.250"	33.438"	6.375"	
104043-00	6	3	2	74.938"	32.125"	79.500"	6.250"	5.250"	33.438"	6.375"	
For 4.3 square foot cages											
104066-00	2	1	2	34.938"	32.125"	79.500"	6.250"	5.250"	33.438"	6.375"	
104070-00	4	2	2	60.968"	32.125"	79.500"	6.250"	5.250"	33.438"	6.375"	
104071-00	6	3	2	86.968"	32.125"	79.500"	6.250"	5.250"	33.438"	6.375"	
For 6 square foot cages											
104122-00	2	1	2	44.938"	32.125"	83.500"	6.250"	5.250"	33.438"	6.375"	
104126-00	4	2	2	80.938"	32.125"	83.500"	6.250"	5.250"	33.438"	6.375"	

Racks with other cage capacities available

* Specify location of downspout; left end or right end.

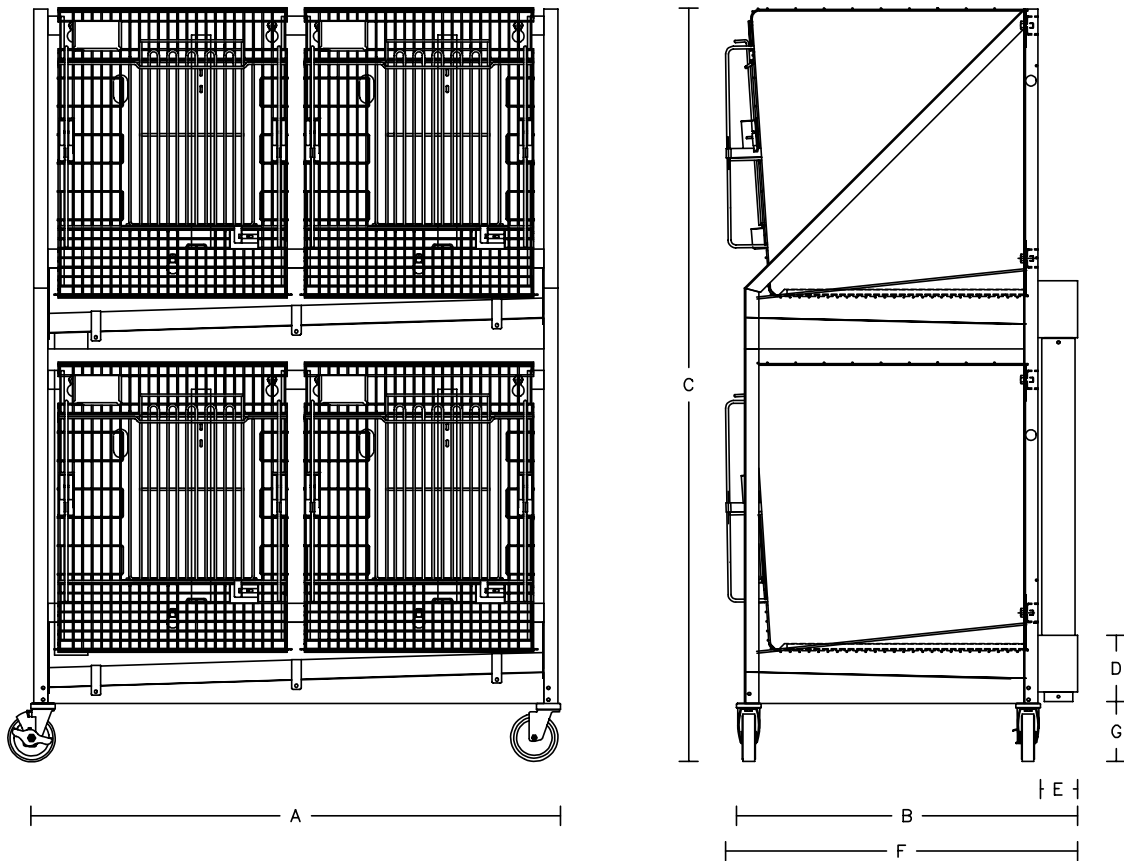

**STAINLESS STEEL ANGULAR FLUSH RACKS WITH BACK DOWNSPOUTS
FOR KEYHOLE SUSPENDED PRIMATE CAGES****

PRODUCT NUMBER	RACK CAPACITY			RACK DIMENSIONS						
	SINGLE WIDTH CAGES	CAGES WIDE	TIERS HIGH	A	B	WITHOUT CAGE C	D	E	WITH CAGE F	G
For 1.6 square foot cages										
104011-00	4	2	2	32.000"	32.625"	66.718"	6.250"	5.250"	33.938"	6.375"
104012-00	6	3	2	48.000"	32.625"	66.718"	6.250"	5.250"	33.938"	6.375"
104013-00	8	4	2	64.000"	32.625"	66.718"	6.250"	5.250"	33.938"	6.375"
For 3 square foot cages										
104031-00	2	1	2	22.000"	39.812"	79.500"	6.250"	5.250"	41.125"	6.375"
104035-00	4	2	2	44.000"	39.812"	79.500"	6.250"	5.250"	41.125"	6.375"
104036-00	6	3	2	66.000"	39.812"	79.500"	6.250"	5.250"	41.125"	6.375"
For 4.3 square foot cages										
104059-00	2	1	2	26.000"	39.812"	79.500"	6.250"	5.250"	41.125"	6.375"
104063-00	4	2	2	52.000"	39.812"	79.500"	6.250"	5.250"	41.125"	6.375"
104064-00	6	3	2	78.000"	39.812"	79.500"	6.250"	5.250"	41.125"	6.375"
For 6 square foot cages										
104115-00	2	1	2	36.000"	39.812"	83.500"	6.250"	5.250"	41.125"	6.375"
104119-00	4	2	2	72.000"	39.812"	83.500"	6.250"	5.250"	41.125"	6.375"

Racks with other cage capacities available

** Specify location of downspout, left rear or right rear.

WALL-MOUNTED SUPPORT SYSTEMS WITH STAINLESS STEEL TROUGHS FOR KEYHOLE SUSPENDED PRIMATE CAGES

Series O Wall-Mounted System supports Series O Cages
Series Q Wall-Mounted System supports Series Q Cages



This Wall-Mounted Support System with Stainless Steel Troughs for Keyhole Suspended Primate Cages, is designed to hang the maximum number of cages on a given length and height wall.

The system is fabricated from electro-galvanized roll-formed channels and trough brackets, with troughs made of stainless steel. Channels are vertically mounted to the wall at both ends and between each cage. Attached to these channels are two horizontal channels per tier of cages. The horizontal channels are fitted with interlocking studs that hold each cage to the system through the keyhole punched slots in the cage backs. Brackets attached to the vertical mounted channels support the stainless steel troughs.

Each trough's ends and back is formed-up above the bottoms of the suspended cages. In addition, the trough's front edge is formed-up to create a lip.

The troughs are custom made to direct waste into a downspout at the center, right end, or left end depending on the drain location in the room.

For additional information on sizing and rack dimensions, please contact SSCI.

These units are built in custom sizes. The dimensions are as built per on-site measurements.

OPTIONS

Stainless Steel watering system
CPVC watering system

STAINLESS STEEL ANGULAR FRONT-TO-BACK FLUSH THROUGH RACKS FOR KEYHOLE SUSPENDED PRIMATE CAGES

Series O Racks support Series O Cages



100005-00 Cages (shown)

104009-00 Rack (shown)

This Stainless Steel Angular Front-to-Back Flush Through Rack for Keyhole Suspended Primate Cages is fabricated from die-formed angles and channels. The angle uprights' lower ends have caster couplers. The angle uprights, angle crossmembers, and channel crossmembers are butted at all joints and welded for strength. Each cage is held securely in place with three keyhole interlocking studs.

Under each tier of cages is a full-width stainless steel trough pitched towards the rear projecting beyond the back of the rack directing waste into various wall troughs or floor gutter systems. The troughs ends are formed up.



The rack has a glass beaded satin finish, removing any discoloration.

The rack is equipped with chrome plated swivel casters with 5" polyolefin wheels, two with wheel brakes. Swivel and axle bearings have zerk grease fittings.

OPTIONS

Stainless Steel watering system
CPVC watering system
Stainless Steel casters
Floor levelers



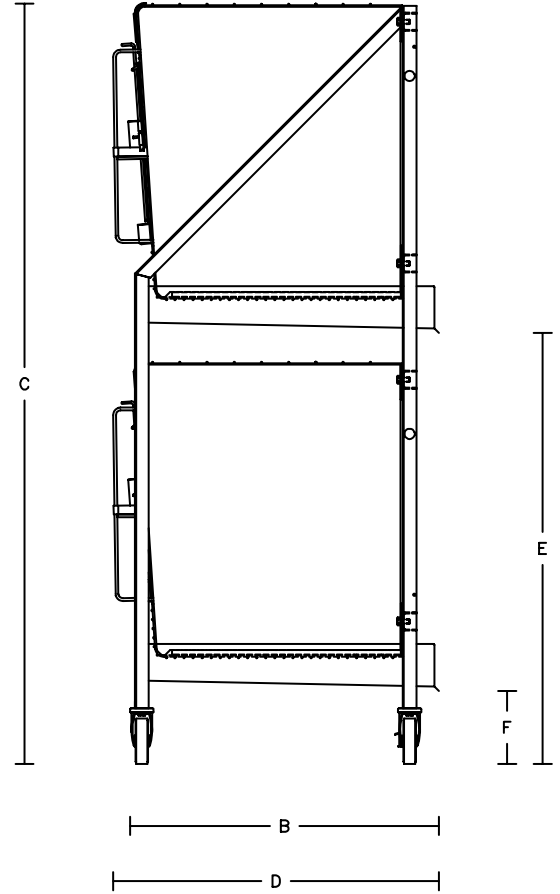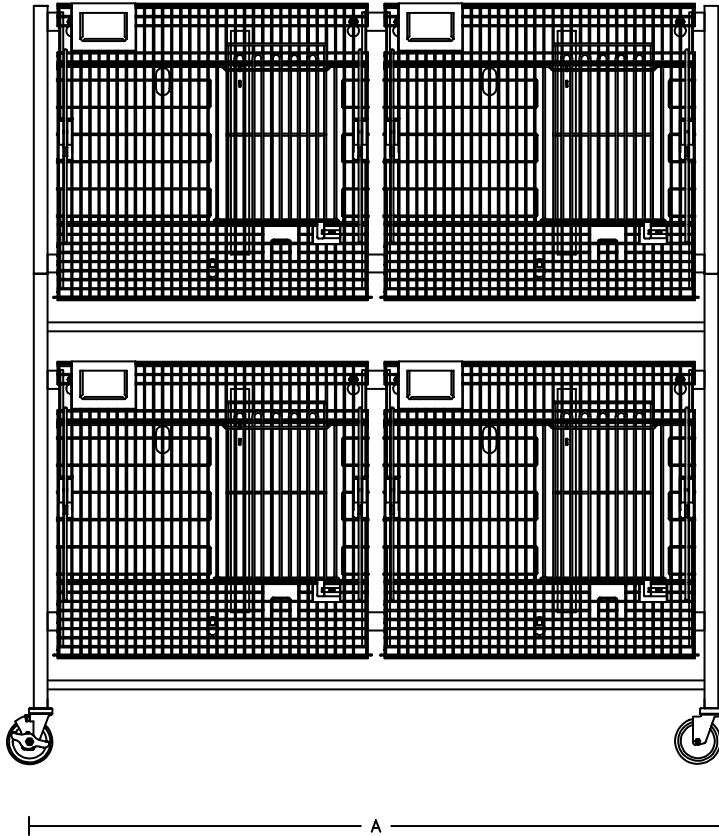

**STAINLESS STEEL ANGULAR FRONT-TO-BACK FLUSH THROUGH RACKS
FOR KEYHOLE SUSPENDED PRIMATE CAGES**

PRODUCT NUMBER	SINGLE WIDTH CAGES	RACK CAPACITY		TIERS HIGH	RACK DIMENSIONS				
		CAGES WIDE	CAGES HIGH		A	WITHOUT CAGE B	C	D	WITH CAGE E
For 1.6 square foot cages									
104007-00	4	2	2	36.938"	26.938"	66.718"	2.000"	26.250"	
104008-00	6	3	2	52.938"	26.938"	66.718"	2.000"	26.250"	
104009-00	8	4	2	68.938"	26.938"	66.718"	2.000"	26.250"	
For 3 square foot cages									
104024-00	2	1	2	26.938"	34.125"	79.500"	2.000"	33.438"	
104028-00	4	2	2	48.938"	34.125"	79.500"	2.000"	33.438"	
104029-00	6	3	2	70.938"	34.125"	79.500"	2.000"	33.438"	
For 4.3 square foot cages									
104052-00	2	1	2	30.938"	34.125"	79.500"	2.000"	33.438"	
104056-00	4	2	2	56.968"	34.125"	79.500"	2.000"	33.438"	
104057-00	6	3	2	82.968"	34.125"	79.500"	2.000"	33.438"	
For 6 square foot cages									
104108-00	2	1	2	44.938"	34.125"	83.500"	2.000"	33.438"	
104112-00	4	2	2	76.938"	34.125"	83.500"	2.000"	33.438"	