



One Company • Many Solutions

## Installation Guide for Digital Temperature Display



This guide provides a visual, step-by-step representation of how to properly install the digital temperature display for the Labrepco Futura and CliniCool Silver Series 1.5 Cu. Ft. Compact Manual Defrost Freezer

Please contact us if you have any additional questions.



One Company • Many Solutions

A close-up photograph showing a person's hand pulling a white, multi-layered door gasket away from the frame of a white appliance door. The gasket is being lifted from the top edge of the door. A red arrow points from a text box to the gasket. The door has a textured surface with several circular indentations. The hand is wearing a black watch with a metal link bracelet.

Pull the door gasket off from the door frame.

Run the green glycol bottle through the gasket.



epCo



Place the display below the "LabRepCo" logo, and connect the probe to the connector.


BY  
CS  
1-800-538-7520  
11-0714 (L)REP/PCD  
Lab, NJ 08809 - USA  
www.ETREKA.com  
1/13 QCRMA0038



Firmly push the gasket back into the door frame.

Place a sticky pad here.



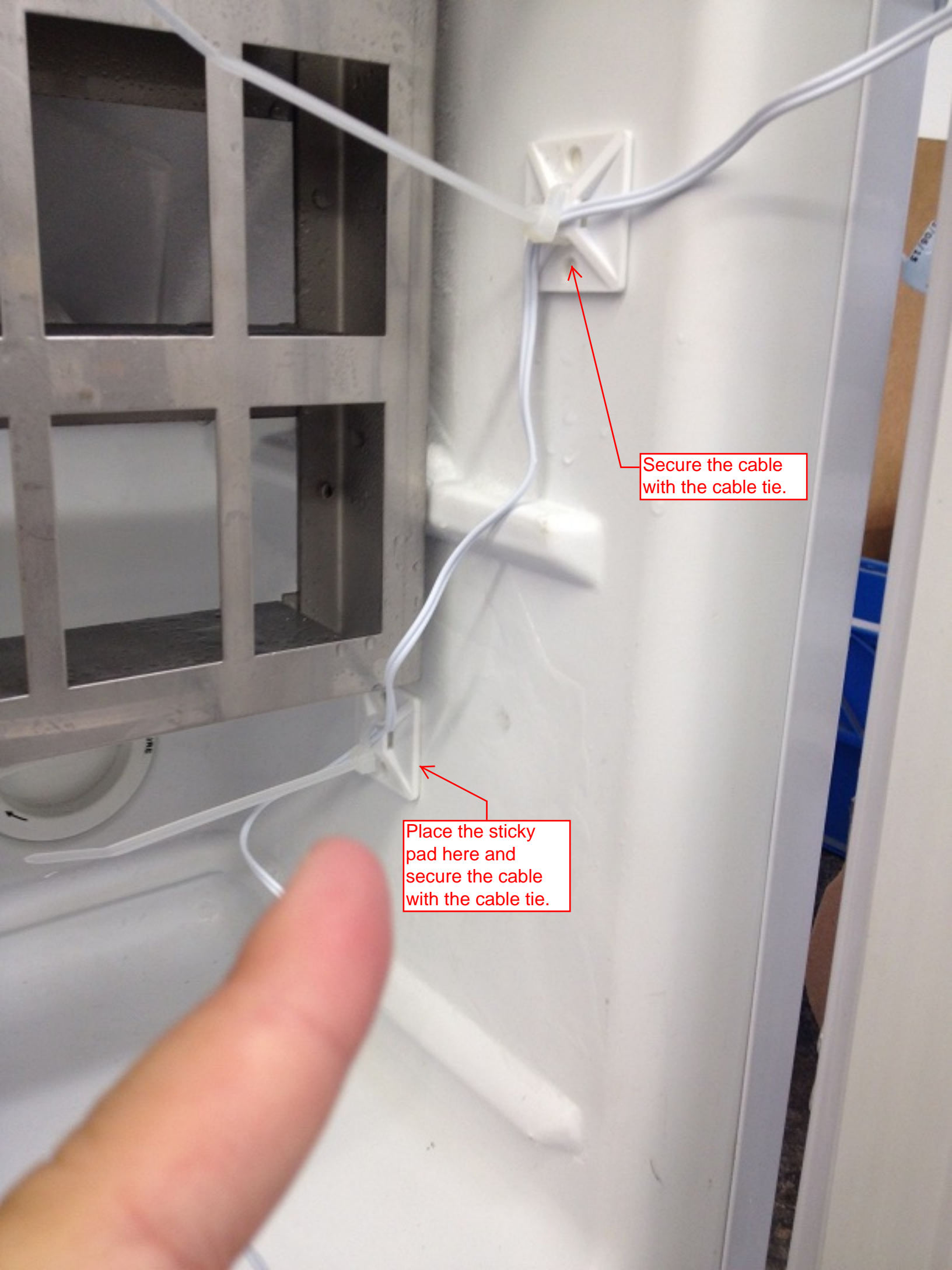

Secure the cable with the cable tie.

Place the sticky pad here and secure the cable with the cable tie.



Roll up the extra cable.






User the cable tie to secure the extra cable.




Remove the backing from the double stick tape on the green glycol bottle.

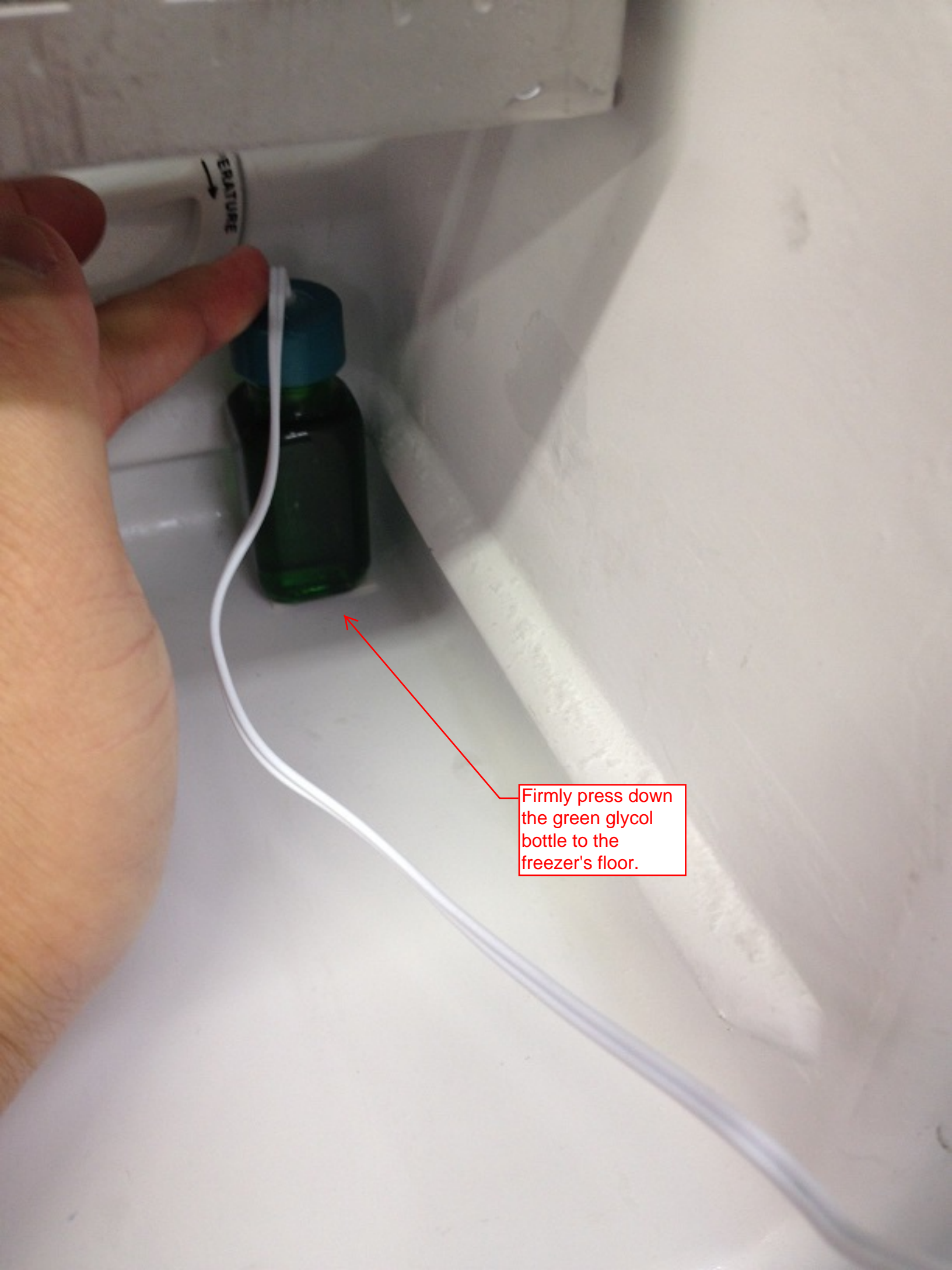

Firmly press down the green glycol bottle to the freezer's floor.

Cut off the extra cable ties off.  
This is the interior completion  
picture.

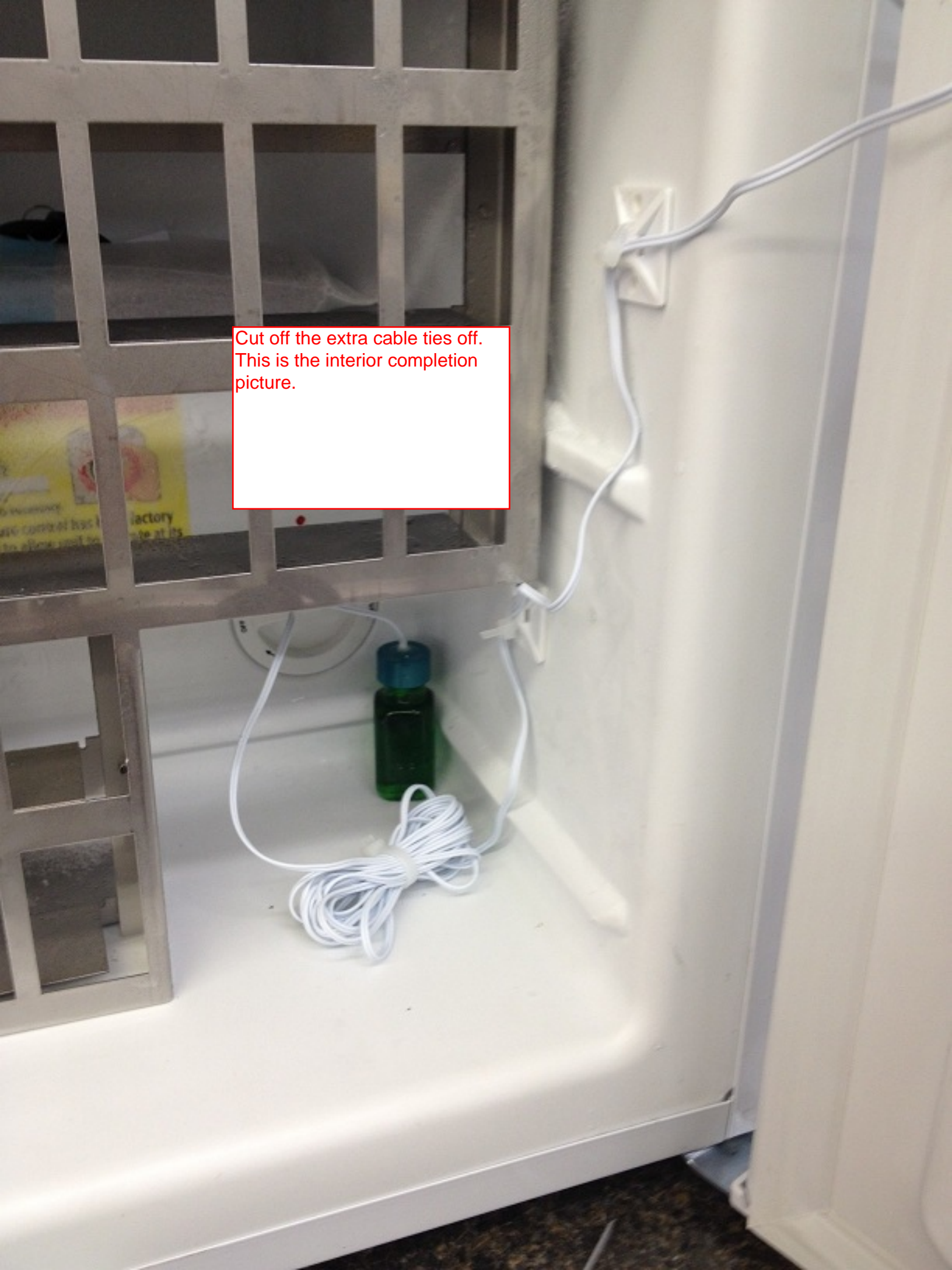

LAB RepCo



This is the exterior completion picture.

800R  
# 130  
LAB REPCO  
# 130  
LAB REPCO