

## **30 cf. shelf reinforcement kit installation**

**HELLO!**

**This kit includes parts to provide additional shelving reinforcement for your 30 cu. ft. freezer.**

**Questions? Feel free to contact us at (800) 648-4041, prompt 5.**

**This kit includes 2 components:**

**1 ea. – Rear Retaining Channel**

**1 ea. – Front Stiffening Channel**

### **STEP 1. Disassembly of the existing shelf**

In order to install the reinforcing components, you must first disassemble the existing shelf by removing the hardware.

1a. Remove and retain 2 existing screws that secure the front of the shelf.



1b. Carefully loosen shelf from existing rear support brackets and remove and retain 2 screws, discard the brackets.



## **STEP 2. Installation of the Rear Retaining Channel**

Using the 2 screws from the existing rear brackets, install the Rear Retaining Channel.



### STEP 3. Installation of the Front Stiffening Channel

Using the screws from front shelf supports, install the Front Stiffening Channel, starting both screws before tightening. Once both screws are partially secure, fully tighten the screws.

