

ULT25NEU Quick Guide:


-----clip, laminate, keep with freezer-----

INITIAL SETUP

Port Stopper/Sealant	!	The Port Stopper/Sealant must be in place during operation.
Power Indicator	A	The unit is ON when lit
Power ON	B	Press and hold for two beeps
Power OFF	B	Press and hold for three beeps
LCD Display	C	Displays chamber temperature
Display Setpoint	F	Press/Release to display current setpoint, S ## displayed
Return to Temp Display	F	Press/Release or wait eight seconds
Change Setpoint	F	Press and hold until <i>Set T LED</i> (E) glows, S ## displayed, then ▲ or ▼ (D)

ADVANCED FUNCTIONS

Change to Preset	Press and hold <i>Set</i> (F) until <i>Set T LED</i> (E) glows, S ## displayed, then select <i>Preset 1, 2 or 3</i> (J)
Postpone Over Temp Alarm	When <i>Over Temp LED</i> (I) glows, press <i>Preset 3</i> (J) until time shows (1.0 h) (If <i>Over Temp LED</i> is not glowing, then <i>Preset 3</i> acts as a Temp Preset)
Return to Temp Display	Wait eight seconds, press <i>Preset 3</i> to return to temperature display
See Error Code Again	If <i>Error LED</i> (H) is glowing, press <i>Preset 2</i> (J) briefly (If <i>Error LED</i> is not glowing, then <i>Preset 2</i> becomes Temp Preset)
Lock Panel	Press and hold <i>Preset 1 AND 2 AND 3</i> (J) until <i>Locked LED</i> (G) turns on (~five seconds)
Unlock Panel	Press and hold <i>Preset 1 AND 2 AND 3</i> (J) until <i>Locked LED</i> (G) turns off (~five seconds)
LCD Display Codes	P ## – Preset Temp, ## h – Alarm Silenced, E ## – Error Code (see operating manual)

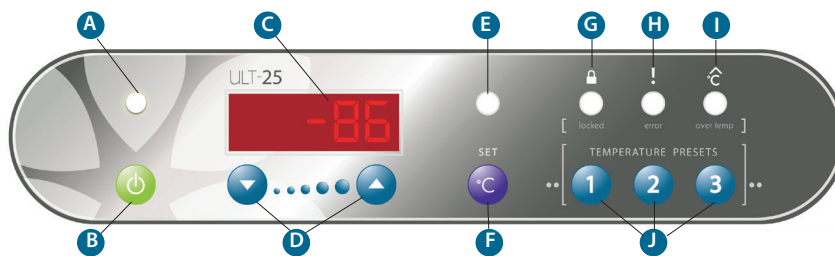

- When unit is first turned on, red Over T light will be on until temperature is within 10° of set point.
- Panel cannot be locked until temperature is within 10° of set point.



6000 Poston Road Athens, Ohio 45701, USA t: 740.274.7900 / 855.274.7900 f: 740.274.7901 www.stirlingultracold.com

©2019 Stirling Ultracold, Global Cooling, Inc. All Rights Reserved.
Global Cooling technology is manufactured under U.S. and International patents.
Stirling Ultracold is a trademark of Global Cooling, Inc.
Specifications subject to change without notice. Refer to www.stirlingultracold.com for the latest specifications.

