

How to Stack a Futura Silver Series Combo Refrigerator/Freezer (L2X Prefix)

In order to properly stack the Combo Refrigerator/Freezers you will need:

- Rubbing Alcohol on a paper towel or soft cloth
- Two pieces of wood or something sturdy enough to hold the weight of the unit
- Stacking kit brackets



Step 1: Clean the surface of the unit that will be on the bottom, with rubbing alcohol



Step 2: Place the two pieces of wood on the surface. Place one centered on the front on the unit and one centered on the back



Step 3: With a partner, lift the unit onto the desired location on top of the pieces of wood.



Step 4: Peel off the adhesive backing located on the stacking kit brackets.



Step 5: Slide the brackets onto all 4 feet of the unit. **Make sure the adhesive is facing down.**



Step 6: Lift the front end of the top unit and remove the piece of wood and carefully place unit down into place. Repeat on the back end of the unit.



Step 7: Check the placement of the unit and firmly push down.

