




Traceable® Thermohygrometer with Clock

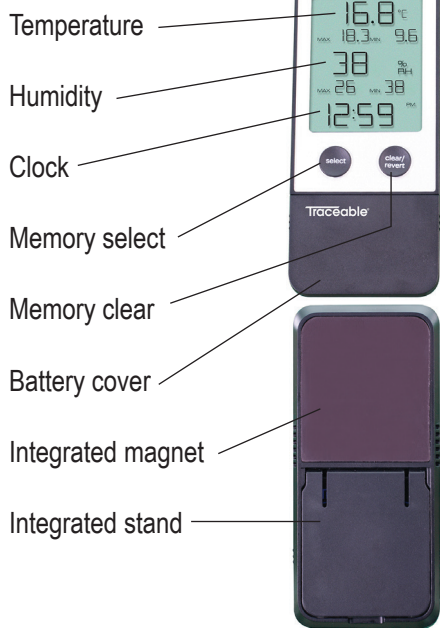

INITIAL SETUP

- 1 Remove battery cover and insert batteries.

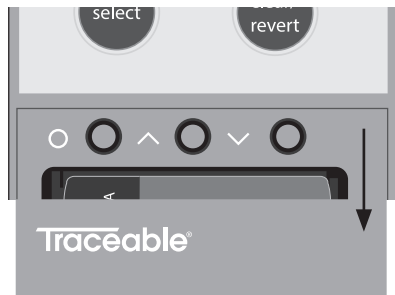

- 2 Once batteries are installed, press  button to select setting.




- 3 The first setting, Hours, will flash on display.



- 4 Press “^” and “v” buttons to adjust setting.




- 5 Once finished, press  button to save setting and proceed to Minutes, 12/24 Hour, etc.




- 6 Complete all settings to configure device:
 - Hours
 - Minutes
 - 12 or 24 hour format clock
 - Temperature units

- 7 Replace battery cover.

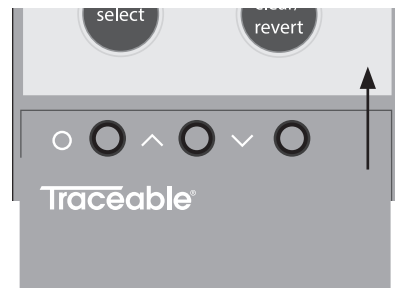

DISPLAY DETAILS

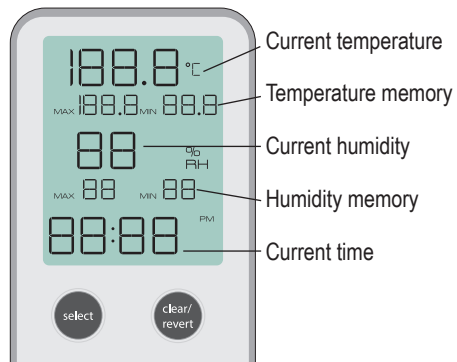

CLEARING INDIVIDUAL READINGS

1. Press select button until desired reading to be cleared is flashing.
2. Press and hold clear/revert button for 3 seconds to clear reading.

CLEARING ALL MEMORY READINGS

1. Press and hold clear/revert button for 3 seconds to clear all memory readings.

WARRANTY, SERVICE, OR RECALIBRATION

Traceable® Products are ISO 9001:2018 Quality-Certified by DNV and ISO/IEC 17025:2017 accredited as a Calibration Laboratory by A2LA.

Traceable® Thermohygrometer with Clock Quick-Start Guide

