TABLEWORX

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⚠ WARNING - TIPPING HAZARD

- ▶ Do not stand, climb, or sit on table.
- ▶ Place heavy objects on lowest shelf.
- ▶ Distribute weight as evenly as possible.
- ▶ Do not drop heavy objects on worktable or subject unit to shock such as forklifts or dollies
- ▶ Move mobile units along long axis (long direction) only (↔).
- ▶ Mobile Units: Bottom structure (solid shelf, three sided frame, H-Frame, etc.) must be on bottom groove.
- ▶ Casters should remain locked except when the cart is in motion.
- ▶ Stationary Units: Bottom structure (solid shelf, three sided frame, H-Frame, etc.) must be as low as possible (NSF requires the lowest shelf be at least 6" from the floor)
- ▶ Riser System: Counterweights **must** be installed for any unit utilizing a Riser System or any overhead accessories.
- ▶ Do not lift by any overhead or bolt-on accessory.
- ▶ Do not lift Table Unit with Work Surface holding from side flanges for shifting purpose

Failure to follow these instructions could result in personal injury or property damage.

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⚠ CAUTION: WEIGHT CAPACITY

650 lbs. (295kg) maximum uniformly distributed static load on the table surface.
 Overhead Accessories: Not to exceed 300 lbs (136kg) total.
 Solid Bottom Shelf: 300 lbs (136kg) maximum uniformly distributed static load per solid bottom shelf.

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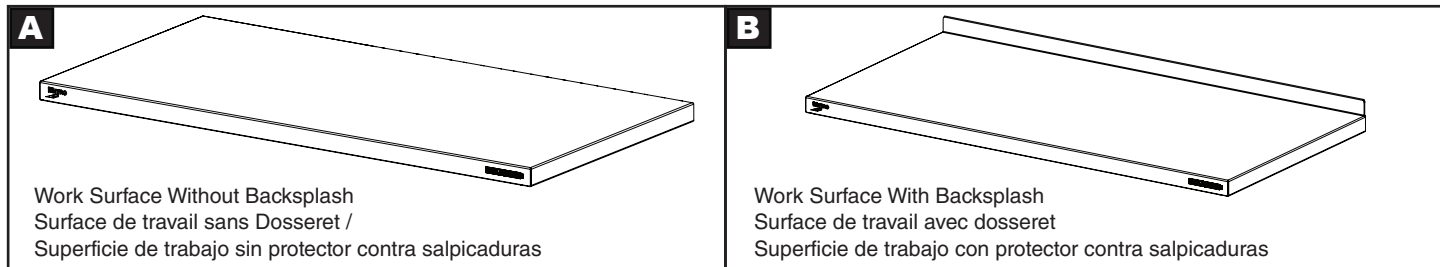
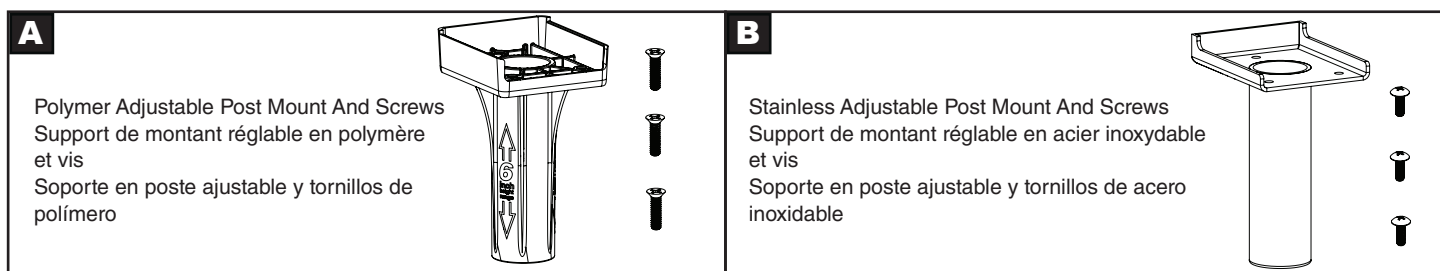
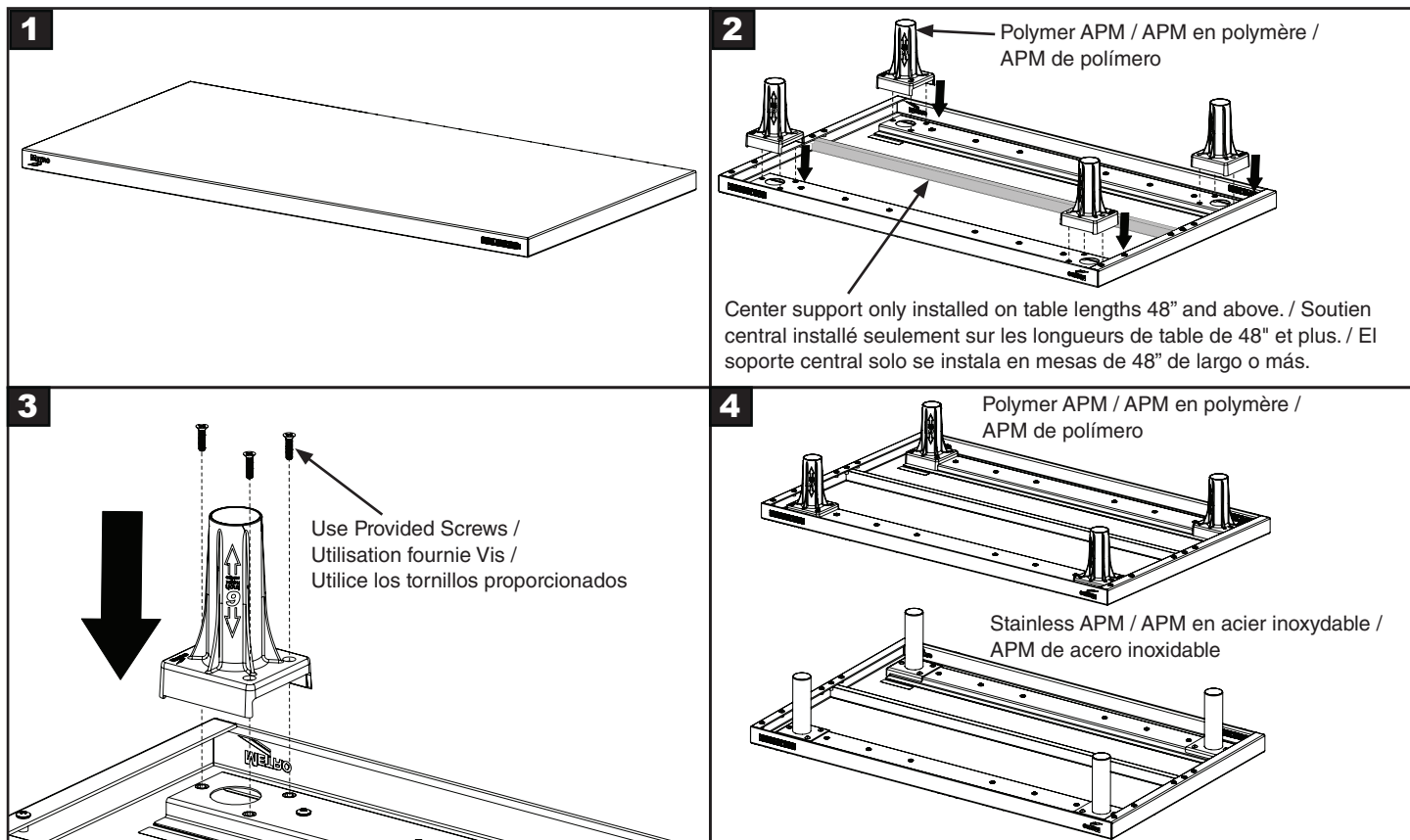
6

⚠ For products with a TableWorx Riser, refer to INSTR-TWRCNTRWT TableWorx Riser for proper installation of **required** counterweights. Scan QR code for instruction.

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4 **⚠ For products with a TableWorx Riser, refer to INSTR-TWRCNTRWT TableWorx Riser for proper installation of required counterweights. Scan QR code for instruction.**

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6

6" (152mm) must be maintained between shelf bottom and floor for proper cleaning

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If necessary, level unit by adjusting leveling feet.

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